

DIMETHYL SULFOXIDE

PHYSICAL PROPERTIES

DMSO



Corporate Office

106 Galeria Blvd.
Slidell, LA 70458
(985) 649-5464

Manufacturing and Technology Support

1880 Fairlawn Rd
Tuscaloosa, AL USA 35401
(205) 561-5023

For samples or further
information related to dimethyl
sulfoxide (DMSO), please contact:

Gaylord Chemical Company, LLC
Customer Service
(800) 426-6620

For more information about
Gaylord Chemical Company, LLC
please visit our website at:
www.GaylordChemical.com

Introduction

Dimethyl Sulfoxide (DMSO) is a highly polar and water miscible organic liquid. It is essentially odorless, and has a low level of toxicity. DMSO is a dipolar aprotic solvent, and has a relatively high boiling point.

This bulletin is a compilation of Physical Properties data for this useful solvent.

Contents

Typical DMSO Properties	page 2
Vapor Pressure vs. Temperature Curve for DMSO	page 3
Specific Gravity of DMSO as a Function of Temperature	page 4
DMSO Viscosity as a Function of Temperature	page 4
Initial Sorption Rates of DMSO at Various Relative Humidities	page 4
Hygroscopicities of DMSO at Various Relative Humidities	page 4
Freezing Temperatures for DMSO - Solvent Mixtures	page 5
Freezing Point Data for DMSO - Water Solutions	page 5
Vapor - Liquid Equilibrium for DMSO - Water Solutions	page 6
Heat Capacity and Density of DMSO	page 7
Boiling Point/Temperature Curves: DMSO - Water Solutions	page 8
Heat of Mixing, DMSO - Water mixtures, 32°C	page 9
Specific Gravity of DMSO - Water Solutions	page 9
Viscosity of DMSO - Water Solutions	page 9
Chemical and Thermal Stability of DMSO	page 9
Comparative Viscosities and Evaporation Times of Common Solvents	page 14



Typical DMSO Properties (in Alphabetical order)

Figure 1

Parameter	Value
Auto ignition temp, in air	300-302°C (572-575°)
Boiling point (1 atmosphere)	189°C (372°F)
Coefficient of expansion	0.00088/°C
Conductivity (Electrical), 20°C	3x10 ⁻⁸ (ohm ⁻¹ cm ⁻¹)
@ 80°C	7x10 ⁻⁸ (ohm ⁻¹ cm ⁻¹)
Critical heat flux (4.10x10 ⁵ J / s / m ²)	1.3x10 ⁵ Btu/hr x ft-2
Critical molar volume	2.38x10 ⁻⁴ m ³
Critical Pressure	56.3 atm. abs.
Critical temperature	447°C (837°F)
Density, at 25°C (see Figure 3)	1.0955 g / cm ³
Dielectric constant, 1 MHz, @ 20°C	48.9
@ 40°C	45.5
Diffusion coefficient	9.0x10 ⁴ cm ² / sec.
Dipole moment, D	4.3
Evaporation rate index @ 25°C	
Relative to n-butyl acetate	0.026
Relative to diethyl ether	0.0005
Flammability limits in air	
lower (100°C)	3 - 3.5% by volume
upper	42 - 63% by volume
Flash point (open cup)	95°C (203°F)
Flash point (closed cup)	89°C (192°F)
Freezing point	18.55°C (65.4°F)
Heat capacity, ideal gas, Cp(T°K)	6.94+5.6x10 ⁻² T
-0.227x10 ⁻⁴ T ²	
Heat capacity (liq.), 25°C	0.47 cal / g / °C
Heat of combustion	6054 cal / g
Heat of fusion	41.3 cal / g
Heat of solution in water at 25°C	-54 cal/g
Heat of vaporization at 70°C	11.3 kcal/mol
Henry's constant @ 21°C	991000
Molar freezing point constant	4.07°C / mol
Molar volume	71.2 cm ³ / g
Molecular weight	78.13
pKa	35.1
pK BH+	-2.7
Refractive index ND@25°C	1.4768

Typical DMSO Properties, continued

Parameter	Value
Solubility parameters	
Hansen's	
- Dispersion	9.0 (cal / cm ³) ^{1/2}
- Polar	8.0 (ca l / cm ³) ^{1/2}
- Hydrogen bonding	5.0 (cal / cm ³) ^{1/2}
Hildebrand's	13.0 (cal / cm ³) ^{1/2}
Specific heat at 29.5°C	0.47 ± 0.015 cal / g / °C
Surface tension at 20°C	43.53 dynes / cm
Vapor pressure at 25°C (See Figure 1)	0.600 mm Hg
Viscosity, cP, at 25°C (See Figure 4)	2.0
Log octanol-water partition coefficient	-1.35

Vapor Pressure vs. Temperature Curve for DMSO

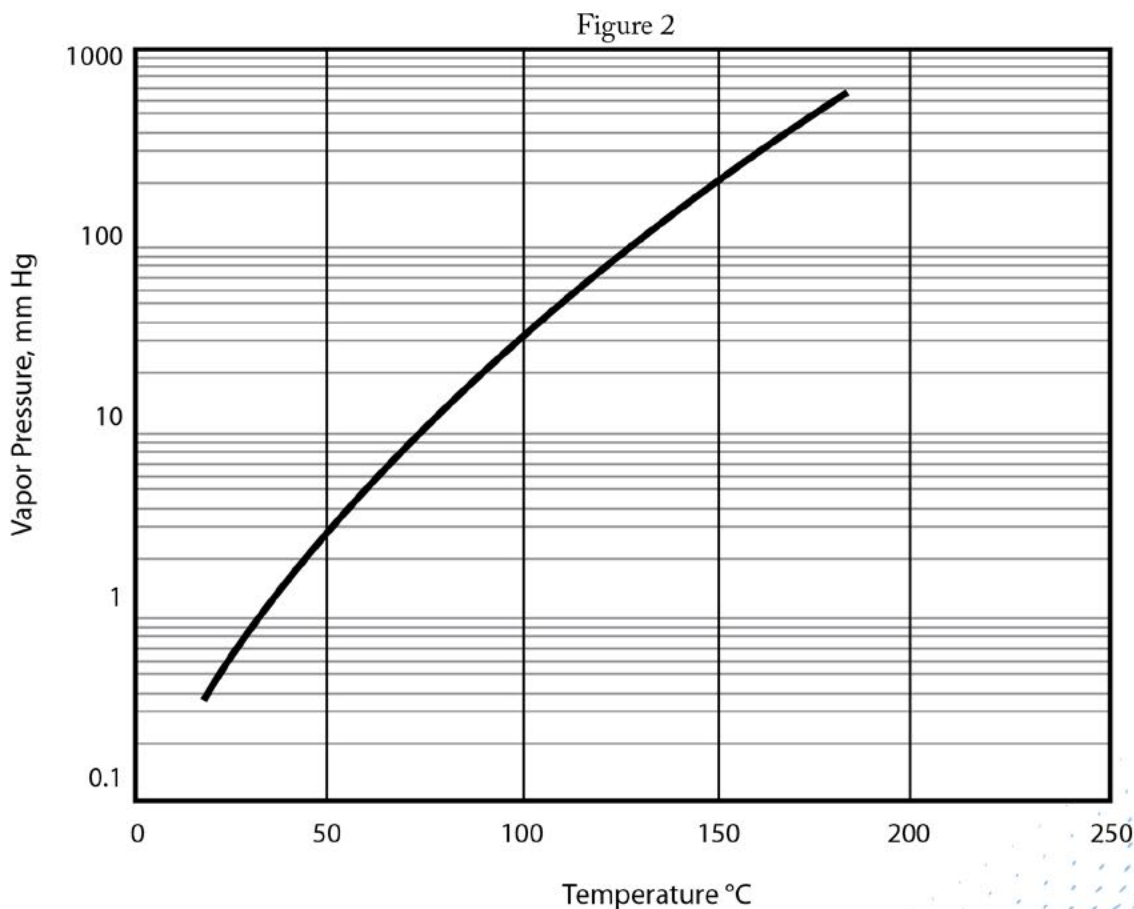




Figure 3

Specific Gravity of DMSO as a Function of Temperature

Temperature (°C)	Specific Gravity (g/cm ³)
15.6	1.1047
21	1.0993
25	1.0955
30	1.0904
40	1.0803
50	1.0702
75	1.0454
100	1.0200
125	0.9946
150	0.974

Figure 4

DMSO Viscosity as a Function of Temperature

Temperature (°C)	Viscosity (cP)
25	1.991
30	1.808
40	1.511
50	1.286
75	0.916
100	0.691
125	0.546

Initial Sorption Rates of DMSO at Various Relative Humidities at 22°C

Figure 5


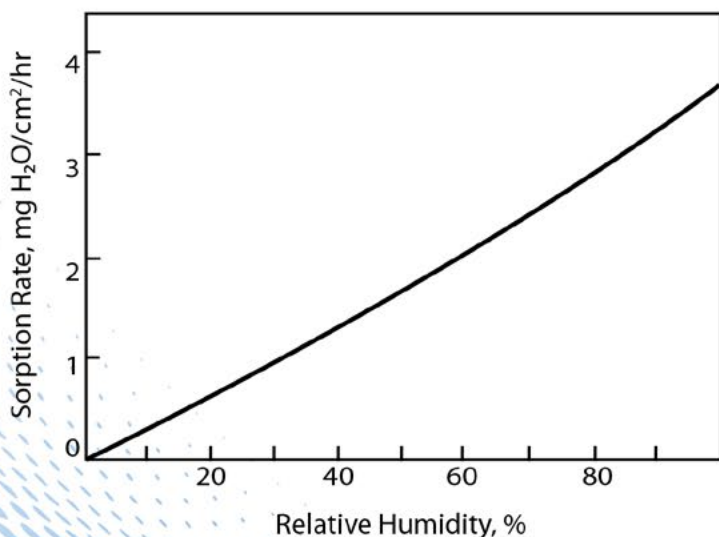
**Hygroscopicities of DMSO at Various Relative Humidities at 22°C**


Figure 6



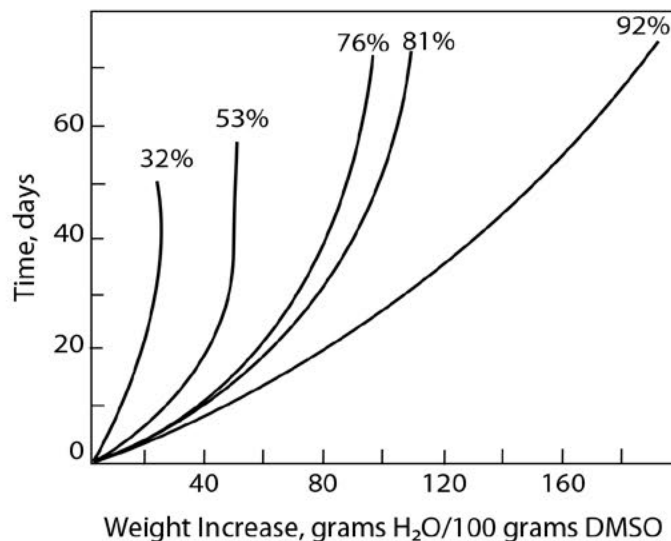

Freezing Temperatures for DMSO-Solvent Mixtures


Figure 7

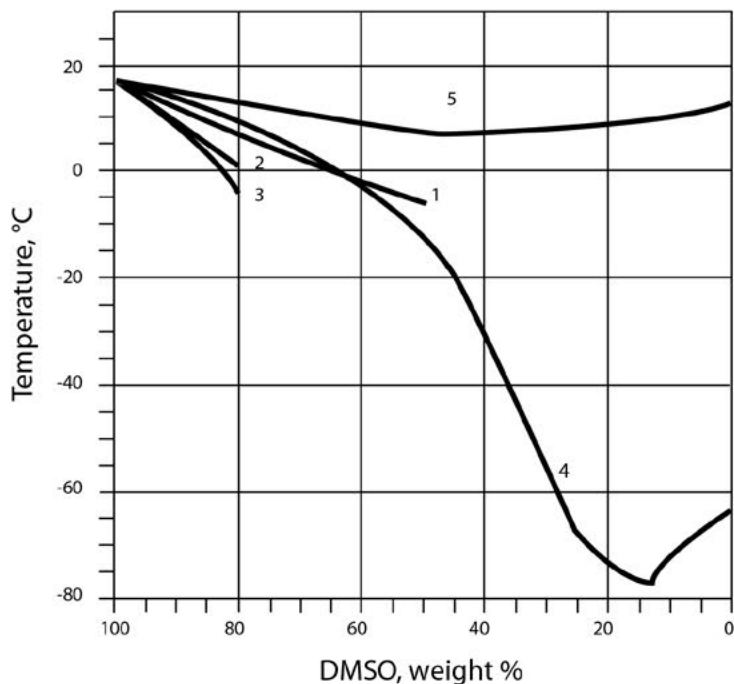

- 1. DMSO - Methyl ethyl ketone
- 2. DMSO - Ethanol
- 3. DMSO - Monoethylene Glycol
- 4. DMSO - Chloroform
- 5. DMSO - p-Xylene

Freezing Point Data for DMSO Water Solutions

Figure 8

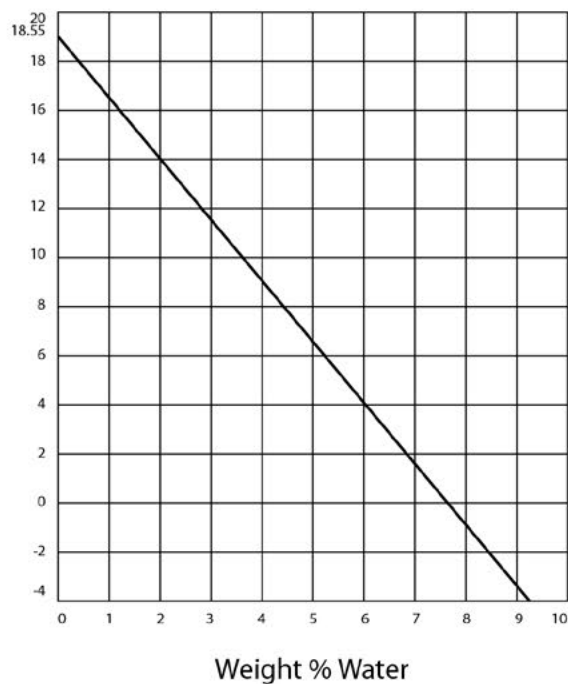

Freezing Point Data for DMSO Water Solutions

Figure 9


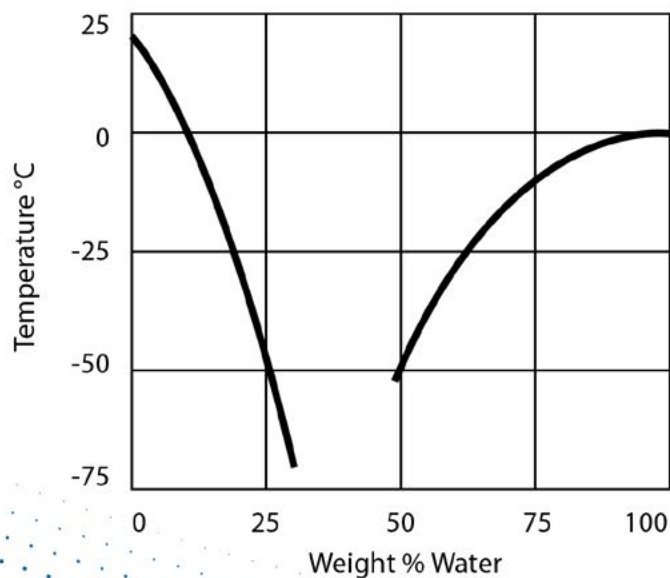


Figure 10



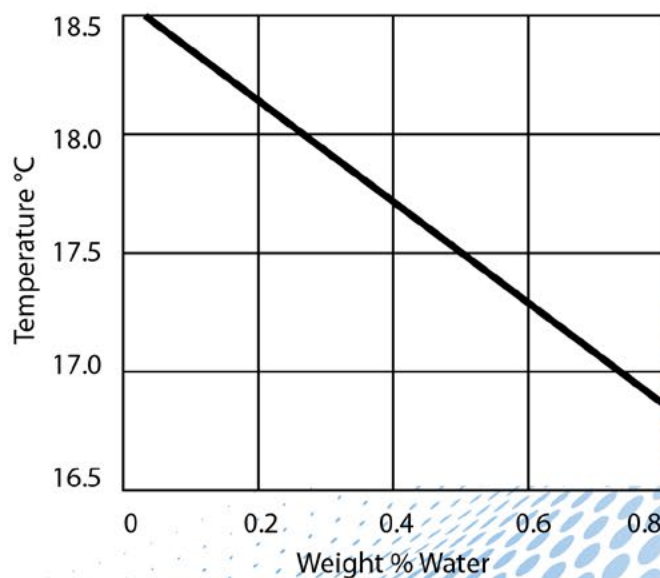

Vapor-liquid Equilibrium for DMSO-Water Solutions (One atmosphere pressure)

Figure 11

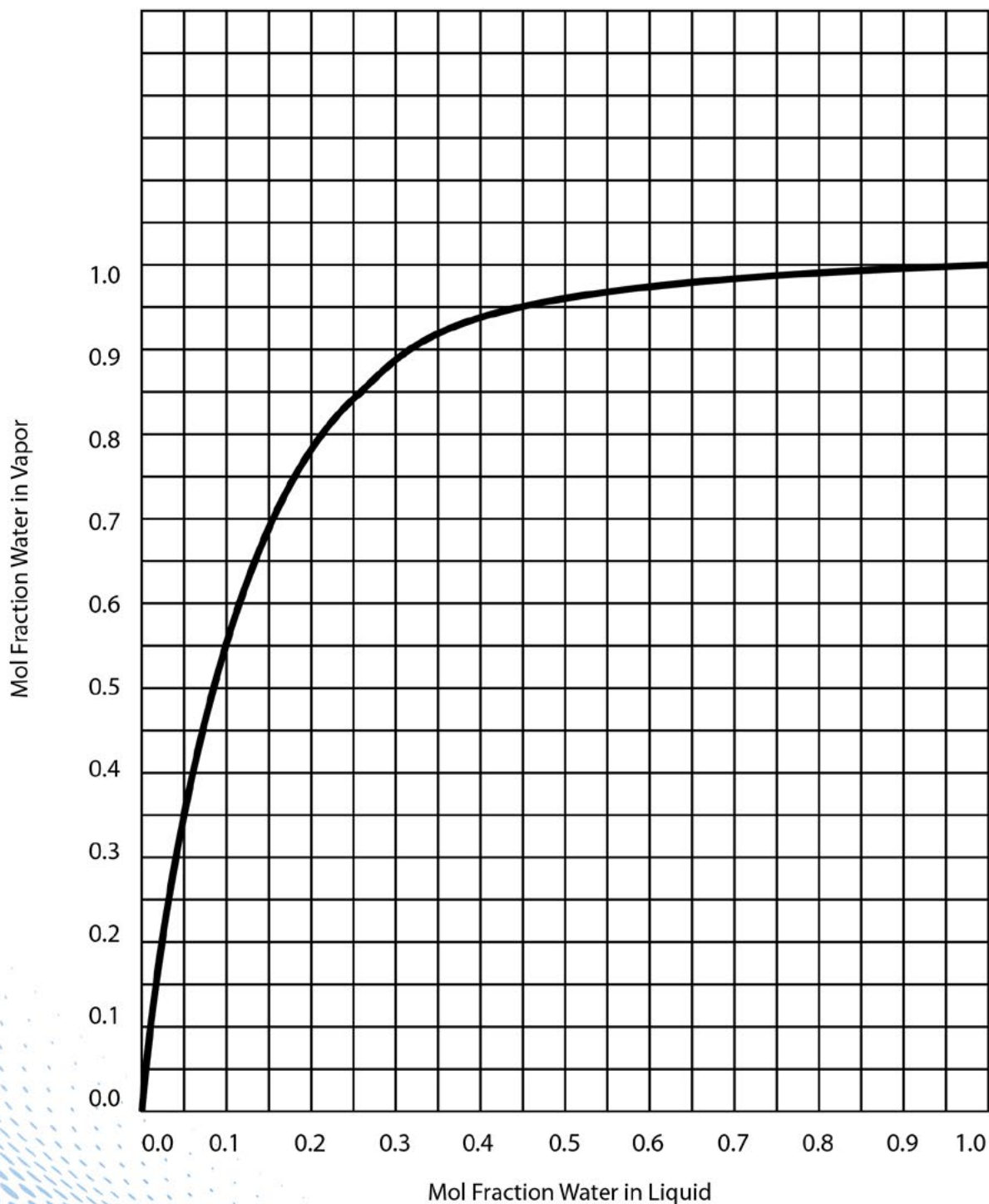




Figure 12

Heat Capacity and Density of DMSO

Temperature, °C	C _p (liquid), cal/(g)(°C)	Density, grams/cc
30	0.47	1.096
60	0.47	1.062
100	0.48	1.023
150	0.52	0.974

Figure 13


**Vapor-Liquid Equilibrium for DMSO-Water Solutions
(One atmosphere pressure)**

Temperature, °C	Mol Fraction Water in Liquid	Mol Fraction Water in Vapor
100.0	1.000	1.000
100.6	0.988	0.9998
101.0	0.975	0.9997
102.0	0.945	0.9994
103.3	0.909	0.9989
105.0	0.865	0.9983
108.0	0.810	0.997
113.0	0.740	0.994
118.0	0.675	0.990
120.0	0.645	0.986
130.0	0.513	0.964
143.0	0.378	0.921
149.0	0.313	0.890
165.0	0.176	0.773
174.5	0.100	0.628
177.0	0.081	0.573
183.0	0.046	0.353
184.6	0.034	0.282
187.7	0.011	0.100
189.0	0.000	0.000




Boiling Point/ Temperature Curves: DMSO Water Solutions (One Atmosphere Pressure)

Figure 14

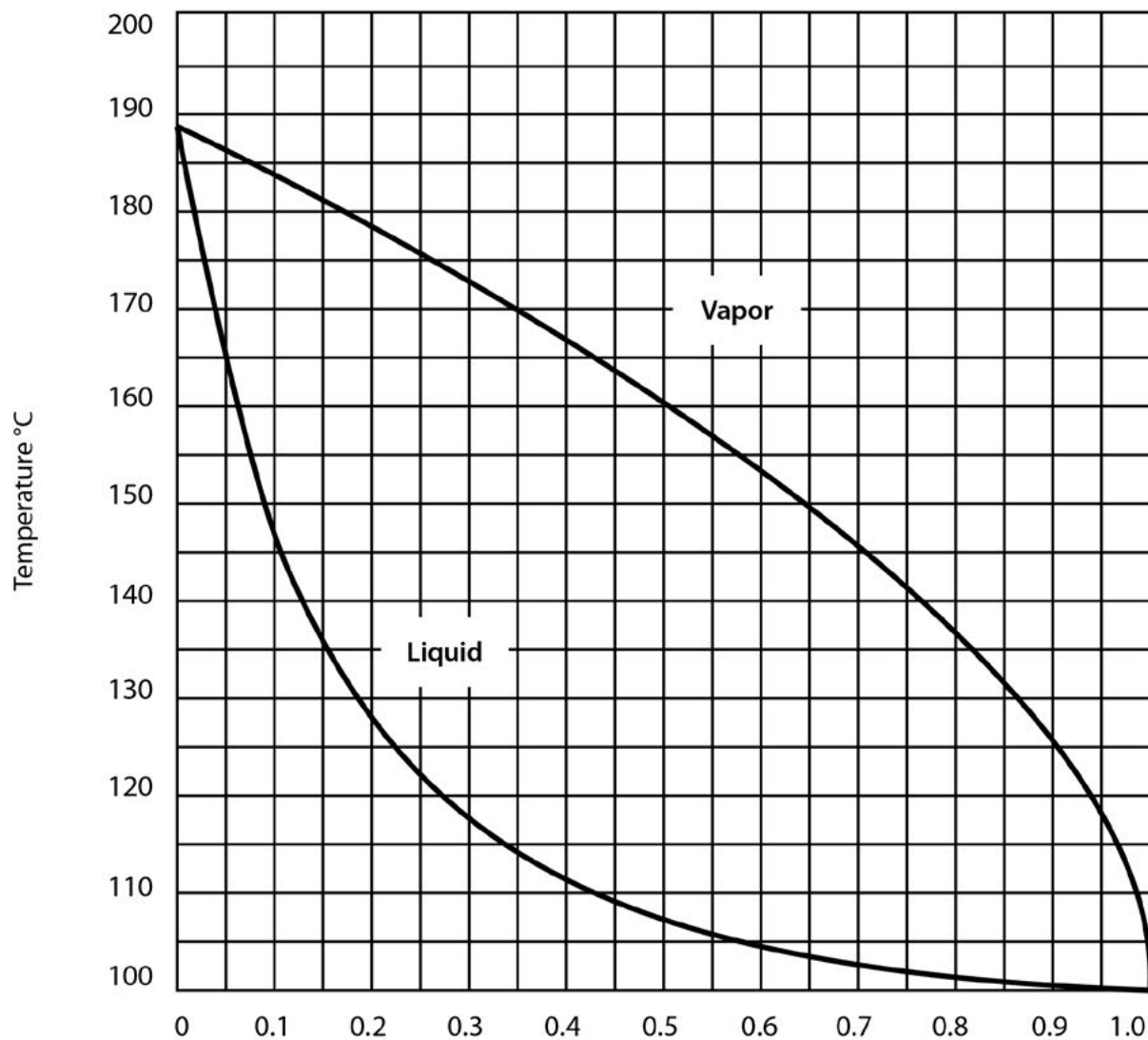


Heat of Mixing of DMSO - H₂O System at 32°C

Figure 15

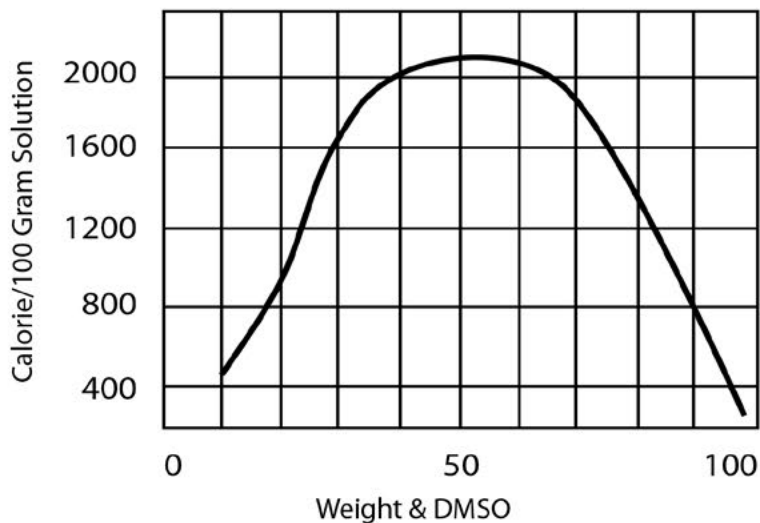


Specific Gravity of DMSO Water Solutions

Figure 16

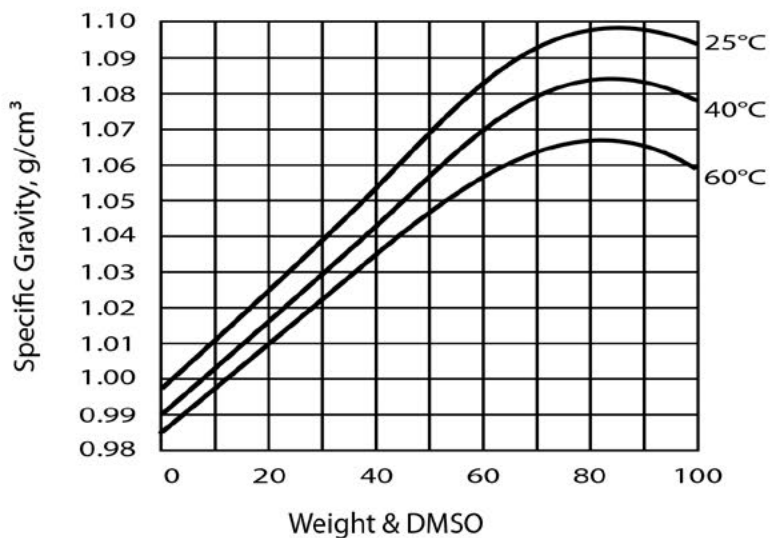

Viscosity of DMSO Water Solutions


Figure 17

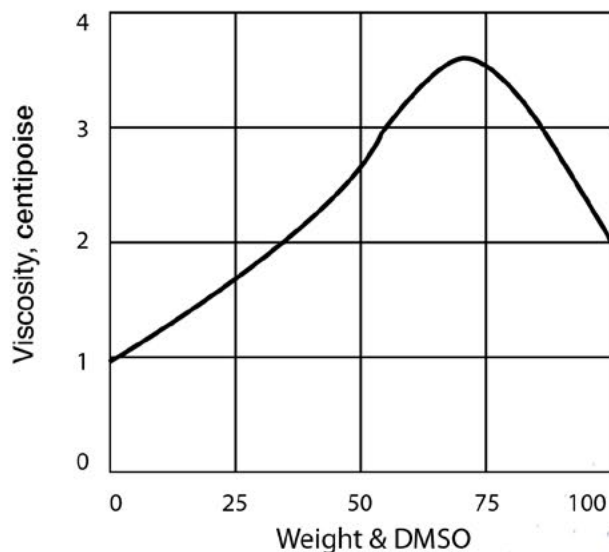

Chemical and Thermal Stability of DMSO

DMSO is highly stable at temperatures below 150°C. For example, holding DMSO at 150°C for 24 hours, one could expect a loss of between 0.1 and 1.0%. Retention times even in batch stills are usually considerably less than this, and therefore, losses would be correspondingly less. It has been reported that only 3.7% of volatile materials are produced during 72 hours at the boiling point (189°C) of DMSO.

Slightly more decomposition, however, can be expected with the industrial grade material. Thus, about 5% DMSO decomposes at reflux after 24 hours. About half of the weight of the volatile materials is paraformaldehyde. Dimethyl sulfide, dimethyl disulfide, bis(methylthio)methane and water are other volatile products. A small amount of dimethyl sulfone can also be found. The following sequence of reactions explains the formation of these decomposition products:

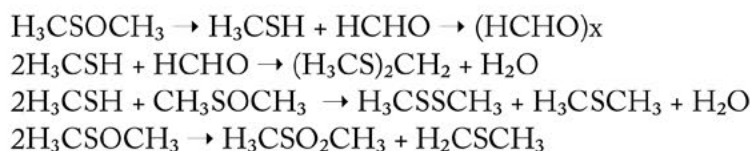

DMSO is remarkably stable in the presence of most neutral or basic salts and bases. When samples of DMSO (300g) are refluxed for 24 hours with 100g each of sodium hydroxide, sodium carbonate, sodium chloride, sodium cyanide, sodium acetate and sodium sulfate, little or no decomposition takes place in most cases. The results are shown in *Figure 18*.

Figure 18
**Results of Reflux with DMSO,
and various materials (24 hours)**

Compound (100g) in 300g DMSO	Reflux Temp., °C	DMSO Recovered % of Original	% Decomposition Products				
			DMS ^(a)	DMDS ^(b)	BMTM ^(c)	HCHO ^(d)	MM ^(e)
NaOH	185-140 ^(f)	93.7	63	31	-	-	-
Na ₂ CO ₃	190	96.3	-	14	-	-	-
NaCl	190	98.7	-	15	-	-	-
NaCN	148-164 ^(g)	100.0	-	-	-	-	-
NaOAc	182-187	97.0	22	33	8	20	-
Na ₂ SO ₄	181-148 ^(h)	85.4	-	-	-	-	-
DMSO Only	189	98.0	15	30	30	-	-

(a) Dimethyl sulfide

(b) Dimethyl Disulfide

(c) Bis-(methylthio)methane

(d) Methyl mercaptan

(e) Formaldehyde


(f) Reflux temperature decreased from 185°C to 140°C over the first 16 hours.

(g) Reflux temperature was 148°C for 20 hours; increased to 164°C during the last 4 hours.

(h) Reflux temperature decreased gradually from 181°C to 148°C.

Thermal Stability of DMSO

Figure 19

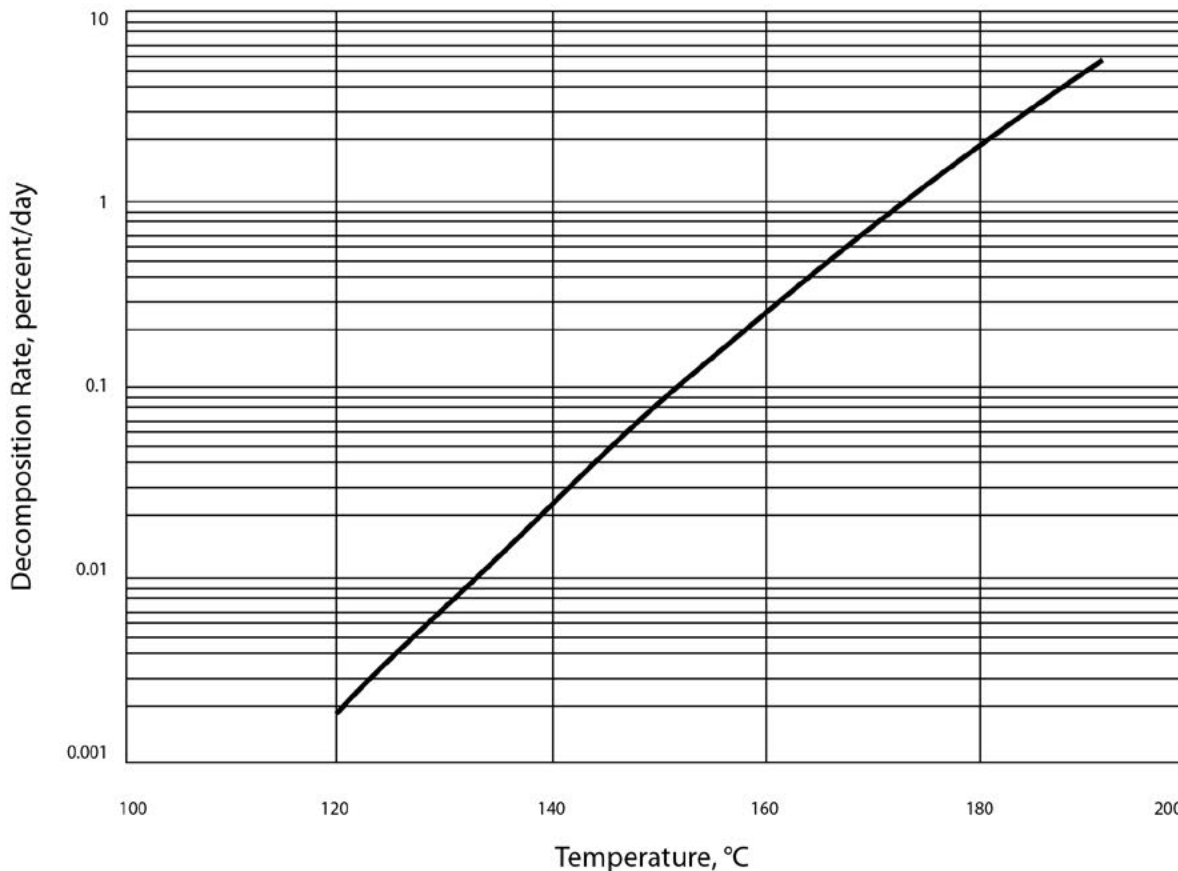

DMSO does not seem to be hydrolyzed by water and very little decomposition of DMSO takes place when it is heated under reflux for periods of 5 to 16 hours. The following tests, shown in Figure 16, have been performed: 1) 10 parts DMSO + 1 part water, 2) 60 parts DMSO + 5 parts water + 1 part sodium hydroxide, 3) 60 parts DMSO + 12 parts water + 1 part sodium bicarbonate, 4) DMSO alone.

Figure 20

Refluxing of DMSO and Mixtures, for shorter periods

Composition of Sample Parts	Reflux Temp., °C	Time, hr	Organic Products Composition, %			
			DMSO	DMS	DMDS	BMTM
10 DMSO: 1 H ₂ O	152	5	100	0	0	0
		15	99.7	0.15	0	0.15
60 DMSO: 5H ₂ O:1 NaOH	155	5	99.8	0.1	0.1	0
		8	99.3	0.6	0.1	0
60 DMSO: 12 H ₂ O: 1 NaHCO ₃	131	6	99.9	0.1	0	0
		12	99.8	0.2	0	0
DMSO Only	191	5	99.8	0.1	0.1	0
		9	99.1	0.2	0.2	0.5
		16	99.0	0.2	0.2	0.6

DMSO is also stable in the presence of concentrated sulfuric or hydrochloric acid at 100°C for up to 120 minutes of heating at atmospheric pressure. Phosphoric acid causes more rapid decomposition of DMSO than does sulfuric or hydrochloric acid. Detected decomposition products are dimethyl sulfide, dimethyl disulfide, and, in smaller quantity, formaldehyde. The results are shown in *Figure 21*:

Figure 21

Effect of heating DMSO with Concentrated Acids

Acid	Conc.	Temp., °C	Time, min.	DMSO % left	% of decomposition product		
					DMS ^(a)	DMDS ^(b)	HCHO ^(c)
H ₂ SO ₄	36N	100	15	99	100		
			30	99	100		
			120	98	100		
H ₂ SO ₄	36N	125	15	86	7	93	
			150	86	7	93	
			210	80	10	90	
H ₃ PO ₄	85%	100	15	92	25	75	
			30	89	45	55	
			45	89	45	55	
			60	87	46	54	
			120	87	46	54	
			150	86	50	50	some
H ₃ PO ₄	85%	125	15	84	25	75	
			60	82	33	67	
			150	82	33	67	

Figure 21

Effect of heating DMSO with Concentrated Acids

Acid	Conc.	Temp., °C	Time, min.	DMSO % left	% of decomposition product		
					DMS ^(a)	DMDS ^(b)	HCHO ^(c)
HCl	12N	95	15	99	100		
			30	99	100		
			60	99	100		
			120	98	100		
HCl	12N	115	15	93	100		
			30	92	100		
			45	87	100		
			60	87	100		
			120	87	100		some

(a) Dimethyl Sulfide (b) Dimethyl Disulfide (c) Formaldehyde

Thermal Decomposition of DMSO

Figure 22

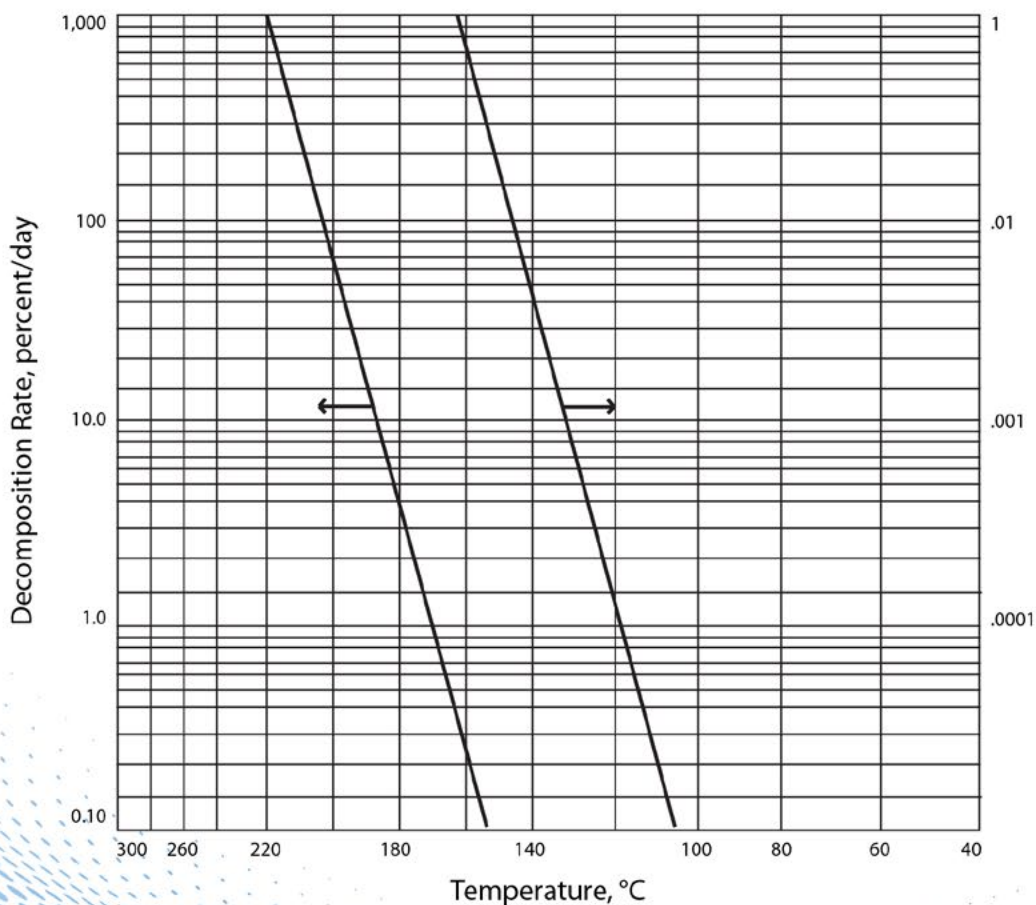




Figure 23

Comparative Viscosities of Common Solvents

Solvent	Viscosity, cps at 25°	
Dimethyl Sulfoxide (DMSO)	2.0	
Dimethylformamide (DMF)	0.8	
N-Methyl-2-pyrrolidone (NMP)	1.6	
Butyrolactone	1.7	
Cyclohexanone	2.1	
Isophorone	2.5	
Diacetone alcohol	3.0	
Propylene Carbonate	4.0	
Sulfolane	10.3	at 30°

Figure 24

Comparative Evaporation Times for Common Organic Solvents

Solvent	90% Evaporation Times, seconds
Dimethyl Sulfoxide (DMSO)	17,600
Dimethylformamide (DMF)	1,570
N-Methyl-2-pyrrolidone (NMP)	2,280
Butyrolactone	3,840
Cyclohexanone	15,400
Isophorone	20,000
Diacetone alcohol	23,700
Propylene Carbonate	119,660
Sulfolane	>1,000,000



**Gaylord
Chemical
Company, L.L.C.**

106 Galeria Blvd., Slidell, LA 70458
(985) 649-5464 (800) 426-6620
www.GaylordChemical.com

The information in this bulletin is based on information available to us and on our observations and experiences. However, no warranty is expressed or implied regarding the accuracy of this data, the results to be obtained from the use thereof, or that any use will not infringe any patent. Each user must establish appropriate procedures for off-loading, handling, and use of the product(s). Since conditions for use are beyond our control, we will make no guarantee of results, and assume no liability for damages incurred by off-loading, handling, or use of the product(s). Nothing herein constitutes permission, or recommendation to practice any invention covered by any patent without license from the owner of the patent.